

## Qualitative tests of Lipids -I-

BCH302 [Practical]

## Lipids

- It can be defined as **nonpolar** organic compound insoluble in polar solvent, but **soluble in organic solvents** such as benzene ,ether, chloroform.
- Lipids are esters of long chain **fatty acids** and alcohols.
- Fatty acids (F.A) are lipids' building blocks.
- There are two types of fatty acids:
- 1. Saturated fatty acids.
- 2. unsaturated fatty acids.

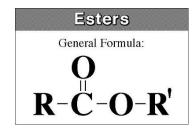


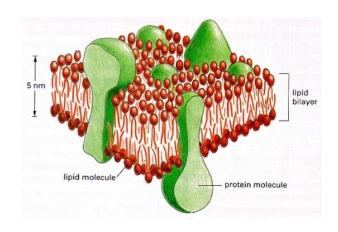


- Biological role of lipids:
- 1. It presents in cell membranes.
- 2. An essential source of energy in the body. It give more energy than carbohydrate and proteins.

## Types of fatty acids



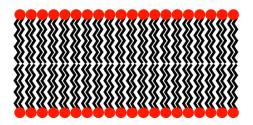

Saturated lipids only

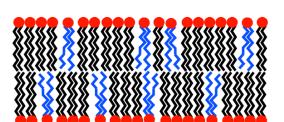


Mixed saturated and unsaturated






Saturated

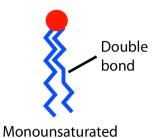

saturated fatty acid

## unsaturated fatty acid

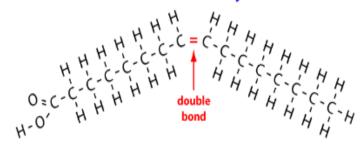

## Classification of Lipids:


- Lipids can be divided according to their chemical composition to:
- I) Simple lipids.
- II) Compound (conjugated) lipids.
- III) Derived lipids.

## I) Simple lipids:

• These compounds are: esters of fatty acids with alcohol.

#### A. Neutral lipids:

Esters of fatty acids with alcohol (glycerol), e.g.: Triacylglyceride.

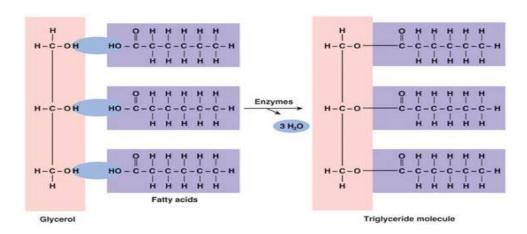

#### **B.** Waxes:

Esters of fatty acids with mono hydroxyle alcohol and higher than glycerol (high molecular weight) e.g.: Beeswax

## II) Compound (conjugated) lipids:

Lipids are linking with <u>other</u> compounds:

## a)Phospholipids:


Most phospholipids contain diglyceride, and phosphate group.


## b)Glycolipids:

Lipids with a carbohydrate attached.

## c)Proteolipids:


Any of a group of proteins to which a lipid molecule is attached.

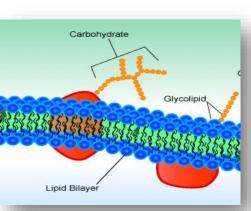

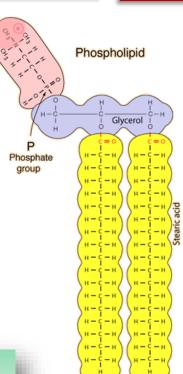

## III) Derived lipids:

• They are substances that are soluble in lipid or derived from the lipids by hydrolysis; for examples, cholesterol and fat soluble vitamins (A, K, E and D).

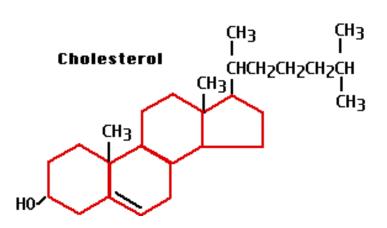



## Practical part

## Qualitative tests of lipids

1 ) Solubility test.

2 Saponification test.

3 Separation of soap from the solution by salting out.

4 Formation of insoluble soap.

## **Experiment 1 : Solubility test**

#### **Objective:**

To test the solubility of oils in different solvent.

## **Principle:**

- Fats are not dissolved in water due to their nature, which is non-polar (hydrophobic), but it is soluble in organic solvents such as chloroform, benzene, and boiling alcohol.
- Different lipids have ability to dissolve in different organic solvent.
- This property enable us to separate a mixture of fat from each other for example:
- ➤ Phosphatidelipid can not dissolve in acetone.
- ➤ Cerebroside and sphingomyline can not dissolve in the ether

## **Experiment 1 : Solubility test**

#### **Method:**

- 1. Place 0.5ml of oil in 6 test tubes clean, dry containing 4ml of different solvents (acetone, chloroform and ether and ethanol, cold ethanol and hot water).
- 2. Shake the tubes thoroughly, then leave the solution for about one minute.
- 3. Note if it separated into two layers, the oil are not dissolve; but if one layer, homogeneous transparent formed, oil be dissolved in the solvent.

Tube	Solvent	Degree of solubility
A	chloroform	
В	ether	
С	ethanol	
D	diluted acid	



dissolved not dissolved

## **Experiment 2 : Saponification test**


• TAG can be hydrolyzed into their component fatty acids and alcohols. This reaction can also be carried out in the laboratory by a process called saponification, where the hydrolysis is carried out in the presence of a strong base (such as NaOH or KOH).

#### **Objective:**

• To form the soap.

#### **Principle:**


- Saponification is a process of hydrolysis of oils or fat with alkaline and result in glycerol and salts of fatty acids (soap).
- Soap can be defined as mineral salts of fatty acids

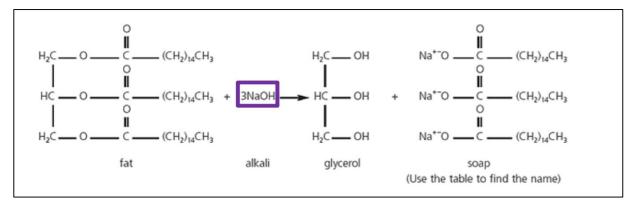



• Soap works on <u>emulsification of oils and fats in the water</u> as it works to reduce the attraction surface of the solution.



• When soap molecules are dissolved in water, the water-repelling hydrophobic tails cluster together while hydrophilic heads surround them arranging themselves in a spherical form toward water molecules.

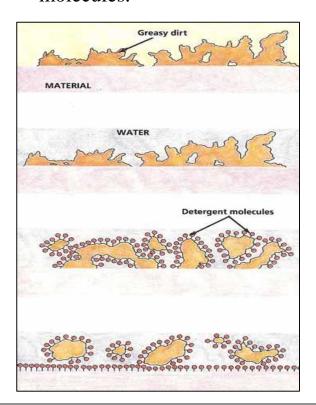

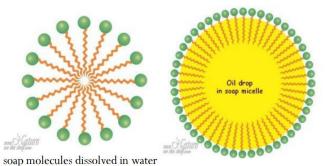


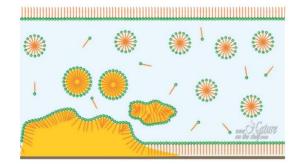

## **Experiment 2 : Saponification test**

#### **Method:**

- 1. Place 2 ml of oil in a large flask.
- 2. Add 4 ml of alcoholic potassium hydroxide.
- 3. Boil the solution for 3 minutes. After this period, make sure it is perfectly saponification process, by taking a drop of the solution and mix with the water if oil separated indicates that the non-completion of the saponification. In this case, continued to boil until all the alcohol evaporates.
- 4. Take the remaining solid material (soap) and add about 30 ml of water and keep it for the following tests.
- 5. Shake the solution after it cools and noted to be thick foam.

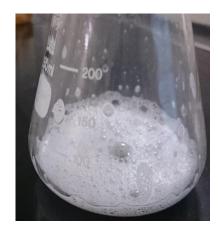

## **Experiment 3 : Testing the separation of soap** from the solution by salting out

## **Objective:**

To investigate the effect of NaCl on soap solubility.

## **Principle:**

- To get the soap out of solution by salting out, when added solid sodium chloride to the solution until saturation; separated soap in the form of insoluble and floats above the surface.
- The NaCl solution provides Na+ and Cl- ions that bind to the polar water molecules, and help separate the water from the soap.
- This process is called salting out the soap.

# **Experiment 3 : Testing the separation of soap** from the solution by salting out

## **Method:**

- 1. Place about 10 ml of soap in the beaker.
- 2. Then add small amounts of sodium chloride in batches, stirring until saturated solution.

Tube	Observation
Soap + NaCl	



## Experiment 4: Test formation insoluble fatty acids salt (insoluble soaps)

## **Objective:**

• To investigate the effect of different cations on soap solubility.

## **Principle:**

- Working calcium, magnesium, lead or iron ions to the **deposition of soap** and make it <u>insoluble in</u> water, where solve these ions **replace the sodium or potassium ions are present in soap**.
- Due to the hard water to contain significant quantities of Ca2+, Mg2+ and some Fe3+ that react with the charged ends of the soaps to form insoluble salts of fatty acid.
- The insoluble salts of fatty acid that Ca2+ and Mg2+ form with soap anions cause white precipitate to from.

 $K^+$  soap +  $Ca^{+2}$  sulfate=>  $Ca^{+2}$  soap +  $K^+$  sulfate.

Potassium soap + calcium sulfate calcium soap + potassium sulfate. (a white precipitate from calcium stearte or oleate is formed).

# **Experiment 4: Test formation insoluble fatty acids salt (insoluble soaps)**

#### **Method:**

- 1. Add about 4 ml of distilled water to 2 ml of soap in two test tubes.
- 2. Add to the first tube a few drops of calcium chloride, to second tube MgCl.

Tube	Observation
CaCl <sub>2</sub>	
MgCl <sub>2</sub>	



## **Howe Work:**

- Why lipids are not soluble in acid?
- What is the chemical composition of your soap?
- Why potassium hydroxide is used in saponification test?